
**Installation instructions
Aufstellanleitung
Notice d'installation
Instrucciones de instalación
Istruzioni di installazione
Installatievoorschrift**

METTLER TOLEDO


**Note
Achtung
Attention
Atención
Attenzione
Opgelet**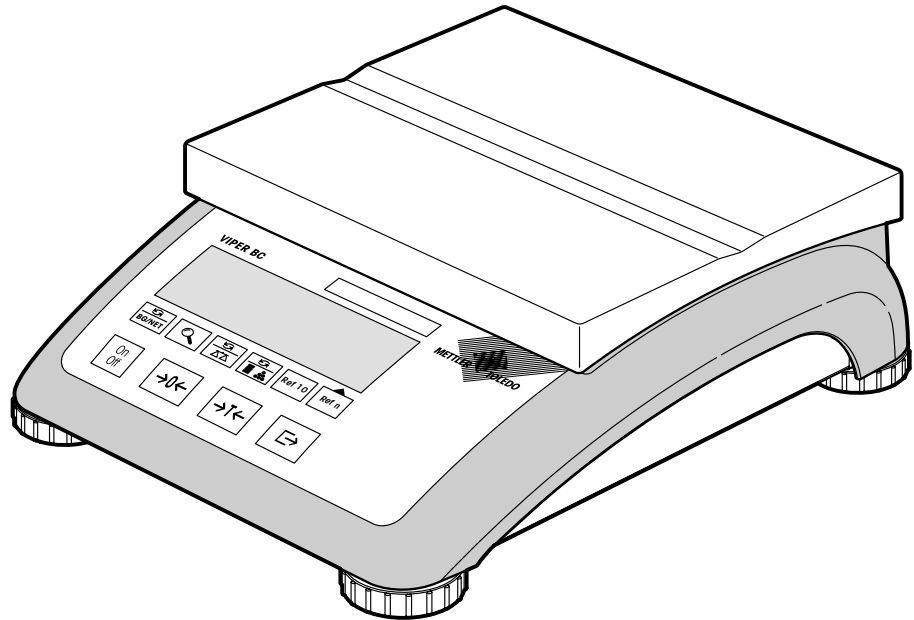

English

Viper scales: The high weighing accuracy, according to the weighing specifications, can only be achieved if the scale is calibrated at its location.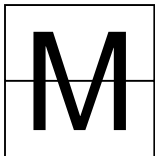

If the packaging bears this mark, the scale or balance has been certified, and all the tests needed to comply with EN45501-8.2 have already been completed at the factory.

Second step of certification at workplace:

The scale or balance must be calibrated at the end user's workplace, and the "metrolo" service mode setting must be changed from non-approved to approved version and the scale or balance switched off and on again.

The green "M" marking must be applied by the responsible weights and measures inspector or authorized service organization.

If national regulations limit the duration of validity of the verification certificate in individual countries, the end user of such a scale or balance is personally responsible for arranging the repeat verification.

For VIPER MonoBloc scales with internal calibration weight (model designation VIPER MBK or VIPER MBDK and designation "IntCal") the following addition to the operating instructions applies:

"IntCal" function: Whenever the power supply is connected, either by plugging the cable into the power supply outlet or (if the scale is battery operated) when the scale is switched on, the scale is automatically correctly calibrated/adjusted using a motorized internal weight. This makes the scale immediately ready for use again if it is moved to a new workplace.

Automatic internal calibration can also be started at any time in master mode using the "CAL" function.

Switch off "IntCal": On non-certified scales in master mode under the second "Scale" block between the "restart" and "Vibrat" menus there is an additional sub-block "ON – CAL" which is set to "ON" at the factory. **The "IntCal" function can be switched off in this sub-block.** (For instructions how to use master mode, refer to the operating instructions, chapter "Scale settings/Scale").

Note: If this setting is set to "Off", the "IntCal" function remains switched off when the scale is connected to the power supply outlet or (if the scale is battery operated) when the scale is switched on.

Special note regarding certifiable or certified scales with "Int-CAL": As a general rule, all scales for EU countries are certified at the factory before shipping. If national regulations in other countries require certification to be done locally, the seal(s) and certification marking(s) are applied by the local weights and measures authorities.

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